

Culturing HEL 299 Cell Line on a Thermo Scientific Nunc Nunclon Cell Culture Treated Surface

Introduction

Thermo Scientific Nunc Nunclon cell culture products are tested for cell growth and plating efficiency using several different cell lines.

Nunc™ Nunclon™ products are tested with two cell lines L929, HEL 299 or F2002 and one Primary Chick Embryo cell culture for monolayer formation, plus cell line V79-4 for cloning efficiency.

The HEL 299 cell line is derived from embryonic lung tissue of a human male. It is a diploid fibroblast-like cell line initially developed for use in vaccine development.

This Tech Note describes a procedure for culturing HEL 299 cell line on a Nunclon treated surface.

Materials and Methods

- HEL 299 cells (ATCC CCL 137)
- Minimum Essential Medium Eagle (MEM)
- Fetal Bovine Serum (FBS)
- L-Glutamine, 200 mM
- Sodium Pyruvate, 100 mM
- Lactalbumin Hydrolysate, 10% in Earle's Balanced Salt Solution
- Dulbecco's Phosphate Buffered Saline, 1X (without Ca²⁺ or Mg²⁺)
- EDTA, 0.02% Solution
- Trypsin Solution, 1X
- Antibiotic/Antimycotic Solution, 100X
- Crystal Violet or Methyl Violet, 0.1-0.4% in aqueous alcohol solution
- Reference lots

Culturing Procedure

1. Place culture vessels (samples and reference lots, as appropriate) in a laminar flow hood along with the culture medium components which have been pre-warmed to 37°C.
2. Prior to harvesting, cells must be at least 75% confluent with good morphology. Aspirate medium and wash cells twice with 1X PBS before addition of EDTA solution.
3. Add EDTA solution to cover the growth area completely.
4. After EDTA solution is decanted, add Trypsin to disaggregate the cells. Incubate culture vessels at 25°C or 37°C, and monitor cell detachment under the microscope. Detachment time will vary.
5. After cells detach, add medium to stop trypsinization and to disperse the cells.
6. Transfer cells to a sterile conical tube and place on ice.
7. Determine cell quantity, e.g. Trypan Blue dye exclusion assay.
8. Determine the number of cells required for each product to be tested by multiplying the plating density by the surface area. Plating density for HEL 299 cell line is 2.0 x 10⁴/cm².
9. Dilute cells into growth medium and seed cell culture product.
10. Incubate cells in a 37°C incubator with 5% CO₂ for seven days to form a confluent monolayer.
11. Decant medium. Add reagent alcohol, 95%, for 5 to 10 minutes for fixation, then decant. Add crystal violet or methyl violet stain, 0.1-0.4%, to cover the surface for 5 to 10 minutes, then decant and wash with water.
12. Evaluate the monolayer when dry (Fig. 1).

Prepare growth medium for HEL 299 cell line as follows:

MEM 1X	500.0 mL
FBS	58.0 mL
L-glutamine	5.8 mL
10% Lactalbumin Hydrolysate	5.8 mL
Sodium pyruvate	5.8 mL
Antibiotic/Antimycotic	5.8 mL
Total	581.2 mL

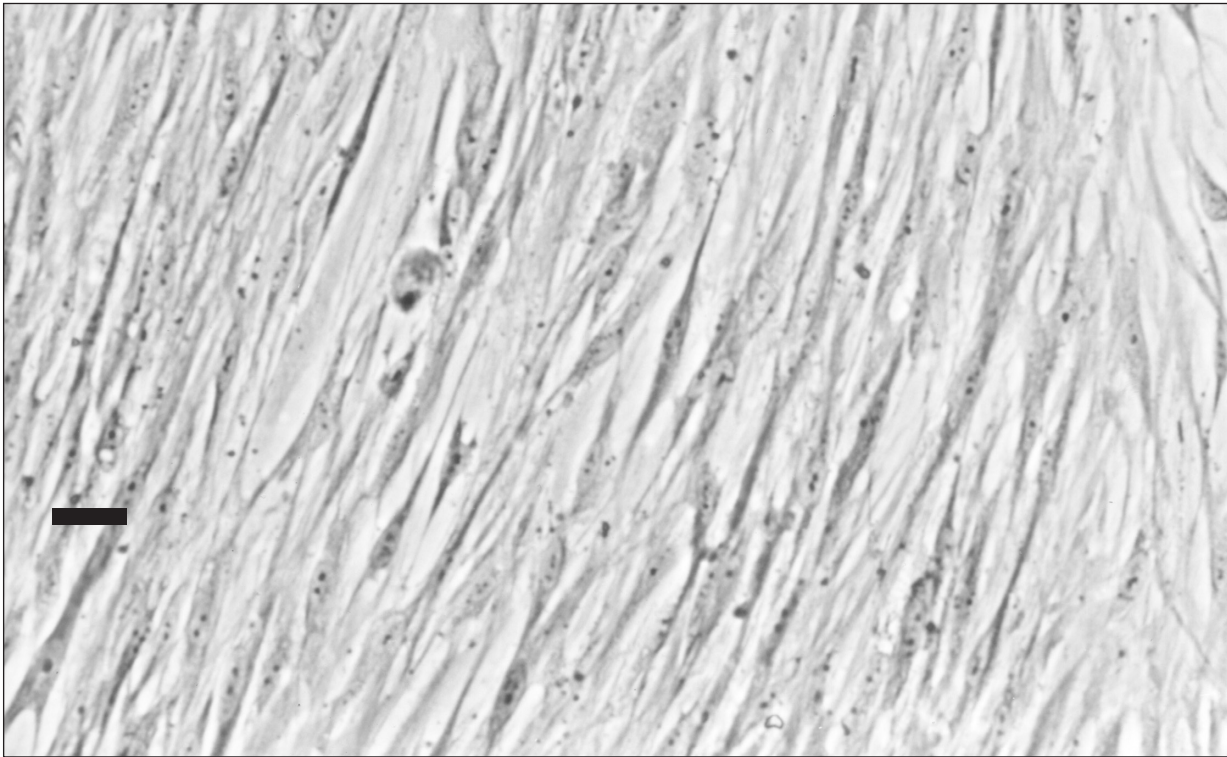


Fig. 1.
HEL299 cells after seven days incubation at 37°C, cultured on a Nunclon polystyrene surface, stained with 0.4% crystal violet. Calibration bar is 40 µm.

Certification Results

When used for Nunclon Certification, HEL 299 cell line results are evaluated as a percentage of surface coverage, per test sample.

- The average percent value must be within 10% of the values of the control products tested with HEL 299 cells.
- Cell growth must be consistent over the entire growth surface.

If these two conditions are met, product passes HEL 299 testing.

Labo Baza

nowoczesne wyposażenie laboratorium

ul. Topolowa 5
62-002 Jelonek k/Poznania
tel.: 061 812 57 45
fax: 061 812 57 25
e-mail: biuro@labobaza.pl
www.labobaza.pl

www.thermoscientific.com

© 2010 Thermo Fisher Scientific Inc. All rights reserved. "Trypsin" is a registered trademark of Sigma-Aldrich Company. All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.

Austria
+43 1 801 40 0

Belgium
+32 53 73 42 41

China
+86 21 68654588

Denmark
+45 4631 2000

France
+33 2 2803 2180

Germany
+49 6184 90 6940

India
+91 22 6716 2200

Italy
+39 02 02 95059 or
434-254-375

Japan
+81 3 3816 3355

Netherlands
+31 76 571 4440

Nordic/Baltic countries
+358 9 329 100

North America
+1 585-586-8800

Russia/CIS
+7 (812) 703 42 15

Spain/Portugal
+34 93 223 09 18

South America
+1 585 899 7298

Switzerland
+41 44 454 12 12

UK/Ireland
+44 870 609 9203

Other Asian countries
+852 2885 4613

Countries not listed
+49 6184 90 6940 or
+33 2 2803 2180

TILSPNUNCTN27 0610